

Ultra Low Temperature Freezer Log

(For Pfizer COVID-19 Vaccine from BAP & VFC programs)

MONTH & YEAR	FREEZER LOCATION/ID	myCAvax ID	VFC PIN

DAY OF MONTH	TIME	INITIALS	ALARM	CURRENT	MIN	MAX	myCAvax/ SHOTS ID
Example	8:00 a.m.	NN		-75.1C	-80.1C	-72.8C	
	4:00 p.m.	NN	✓	-65.2C	-78.3C	-58.8C	12345
1	a.m.						
	p.m.						
2	a.m.						
	p.m.						
3	a.m.						
	p.m.						
4	a.m.						
	p.m.						
5	a.m.						
	p.m.						
6	a.m.						
	p.m.						
7	a.m.						
	p.m.						
8	a.m.						
	p.m.						
9	a.m.						
	p.m.						
10	a.m.						
	p.m.						
11	a.m.						
	p.m.						
12	a.m.						
	p.m.						
13	a.m.						
	p.m.						
14	a.m.						
	p.m.						
15	a.m.						
	p.m.						

Notes: _____


Keep all temperature logs and data files for three years.

VFC Program: (877) 243-8832

BAP: (833) 502-1245

Instructions

Keep ULT freezer in OK range.

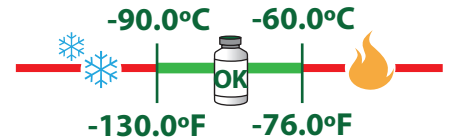

Check temperatures twice a day.

1. Fill out month, year, refrigerator ID, myCAvax ID and VFC PIN.
2. Record the time and your initials.
3. Record a check if an alarm went off.
4. Record Current, MIN, and MAX.
5. Clear MIN/MAX.
6. Ensure data logger is recording.

IF ALARM WENT OFF:

1. Post "Do Not Use Vaccines" sign.
2. Alert your supervisor.
3. Report excursion to myCAvax for BAP vaccine and MyVFCvaccines for VFC vaccine.
4. Record incident ID from myCAvax (BAP) and SHOTS (VFC).

Supervisor's Review

When log is complete, check all that apply:

- Month/year/location/ID/PIN are recorded.
- Temperatures were recorded twice daily.
- I reviewed data files for all the days on this log to find any missed excursions.
Date downloaded: _____
- Any excursions were reported to myCAvax (BAP) or SHOTS (VFC).
- We understand that falsifying this log is grounds for vaccine replacement and termination from the BAP/VFC programs.

On-Site Supervisor's Name: _____

Signature: _____

Date: _____

Staff Names and Initials: _____

Ultra Low Temperature Freezer Log

(For Pfizer COVID-19 Vaccine from BAP & VFC programs)

MONTH & YEAR	FREEZER LOCATION/ID	myCAvax ID	VFC PIN

DAY OF MONTH	TIME	INITIALS	ALARM	CURRENT	MIN	MAX	myCAvax/ SHOTS ID
16	a.m.	-----					
	p.m.	-----					
17	a.m.	-----					
	p.m.	-----					
18	a.m.	-----					
	p.m.	-----					
19	a.m.	-----					
	p.m.	-----					
20	a.m.	-----					
	p.m.	-----					
21	a.m.	-----					
	p.m.	-----					
22	a.m.	-----					
	p.m.	-----					
23	a.m.	-----					
	p.m.	-----					
24	a.m.	-----					
	p.m.	-----					
25	a.m.	-----					
	p.m.	-----					
26	a.m.	-----					
	p.m.	-----					
27	a.m.	-----					
	p.m.	-----					
28	a.m.	-----					
	p.m.	-----					
29	a.m.	-----					
	p.m.	-----					
30	a.m.	-----					
	p.m.	-----					
31	a.m.	-----					
	p.m.	-----					

Notes: _____


Keep all temperature logs and data files for three years.

VFC Program: (877) 243-8832

BAP: (833) 502-1245

Instructions

Keep ULT freezer in OK range.

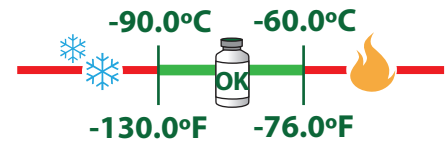

Check temperatures twice a day.

1. Fill out month, year, refrigerator ID, myCAvax ID and VFC PIN.
2. Record the time and your initials.
3. Record a check if an alarm went off.
4. Record Current, MIN, and MAX.
5. Clear MIN/MAX.
6. Ensure data logger is recording.

IF ALARM WENT OFF:

1. Post "Do Not Use Vaccines" sign.
2. Alert your supervisor.
3. Report excursion to myCAvax for BAP vaccine and MyVFCvaccines for VFC vaccine.
4. Record incident ID from myCAvax (BAP) and SHOTS (VFC).

Supervisor's Review

When log is complete, check all that apply:

- Month/year/location/ID/PIN are recorded.
- Temperatures were recorded twice daily.
- I reviewed data files for all the days on this log to find any missed excursions.
Date downloaded: _____
- Any excursions were reported to myCAvax (BAP) or SHOTS (VFC).
- We understand that falsifying this log is grounds for vaccine replacement and termination from the BAP/VFC programs.

On-Site Supervisor's Name: _____

Signature: _____

Date: _____

Staff Names and Initials:

